

'CUSTOM MACROS' PLUGIN INSTALLATION (non-network)

- **ImageJ Site:**
 - <https://imagej.nih.gov/ij/download.html>
 - **'Custom Macros' Plugin Site:**
 - <https://ijmacros.com>
 - This tutorial shows installation in Linux (*Mint 18*).
However, since ImageJ is a portable, cross-platform, Java program, the procedure should be very similar for any operating system.
-

1. Download the ImageJ archive zip file.
2. Extract the ImageJ program to a directory of your choice.
 - (*We will use the desktop for convenience.*)
3. Download the 'Custom Macros' archive zip file.
4. Extract the plugin to /ImageJ/macros/ folder.
 - Overwrite when prompted.
5. Optionally (preferably) create a desktop link to ImageJ.
 - Install recommended plugins (EDF and Stitching).
 - (*Not shown*).
6. Run ImageJ and press the **orange power-on 'Start button'**.
7. Follow instructions on calibration.
 - An example using a camera manual is given.
8. **INSTALLATION IS COMPLETE!**
 - *Network setup is simple and shown in another video.*
9. Please watch the **User Sign-In Tutorial** next.